

Lovibond® Colour Measurement Tintometer® Group



AF330 EBC Colour Scale

2-Field Visual Measurement of Beer According to the EBC Scale



- Measure Beer According to EBC Colour Over the Range 2 - 27
- Fast, Economical & Portable Visual Matching
- Two-Field Vision for Accurate Results
- Ability to Guarantee Constant Lighting Conditions

Part Number: 433300

Ensures Your Beer Production is Always Optimal

By measuring the colour of your beer, you can rest assured that your customers are always receiving the same colour of the drinks, reassuring them of the quality of your produce. The Lovibond® AF 330 is a special kit based on the Lovibond® Comparator 2000+ for visual colour grading according to EBC Colour over the range 2 - 27. The colour is determined by matching the sample in a 25 mm optical glass cell with the precalibrated coloured glass filters.

Fast, Economical & Portable Visual Matching

The Comparator 2000+ is a flexible, modular system for Visual Colour Grading. Using a suitable comparator instrument and glass disc, you can quickly visually match the sample directly against the calibrated, colour stable glass standards.

Two-Field Vision for Accurate Results

Lovibond® comparator instruments include an integrated prism which causes the sample and comparison fields to overlap, allowing enhanced colour discrimination and matching. The prism is ultrasonically sealed to prevent contamination of the optics. Each individually graded glass standard represents a discrete step on the measurement range (as opposed to a continuous graduation), aiding colour discrimination and allowing great accuracy of results.

Ability to Guarantee Constant Lighting Conditions

An optional item for the Comparator 2000+ is the Daylight 2000 Lighting Unit. This is a standardised light source which guarantees you a constant lighting condition for accurate colour grading. This is particularly useful when the sample is very pale in colour. By using a Lovibond® lighting unit, you can be confident of uniform lighting conditions for colour matching, 24 hours a day and irrespective of ambient lighting.

Extended Shelf Life

The unique feature of Lovibond® test discs is that they incorporate coloured glass standards that are colour stable and not affected by UV light or extreme environmental conditions.

Please note, we cannot guarantee compatibility of discontinued Lovibond® Comparator 1000 disks with the Lovibond® 2000+ Comparator.

Industry

Food and Beverage Industry

Application

Beer, Malts and Caramels

AF330 EBC Colour Scale

The Lovibond® AF 330 is a special kit based on the Lovibond® Comparator 2000+ for visual colour grading according to EBC Colour over the range 2 - 27. Grading to a colour scales ensures that your customers always receive the product they expect, inspiring confidence in the quality of the produce.

Technical Data

Applicable Standard EBC Analytika

Colour Type Transparent

Portability Portable

Delivery Scope

- 1 x Lovibond 2000+ Comparator
- 1 x Daylight 2000 Unit (EBC)
- 1 x 4/14A EBC Colour Scale Disc
- 1 x 4/14B EBC Colour Scale Disc
- 1 x 4/14C EBC Colour Scale Disc
- 1 x 4/14D EBC Colour Scale Disc
- 2 x W680.0G.25mm Cells

Tintometer GmbH

Lovibond® Water Testing
Schleefstraße 8-12
44287 Dortmund
Tel.: +49 (0)231/94510-0
Fax: +49 (0)231/94510-30
sales@lovibond.com
www.lovibond.com
Germany

Tintometer China

Room 1001, China Life Tower
16 Chaoyangmenwai Avenue,
Beijing, 100020
Tel.: +86 10 85251111 App. 330
Fax: +86 10 85251001
chinaoffice@tintometer.com
www.lovibond.com
China

Tintometer Inc.

6456 Parkland Drive
Sarasota, FL 34243
Tel: 941.756.6410
Fax: 941.727.9654
sales@lovibond.us
www.lovibond.com
USA

The Tintometer Limited

Lovibond House
Sun Rise Way
Amesbury, SP4 7GR
Tel.: +44 (0)1980 664800
Fax: +44 (0)1980 625412
sales@lovibond.uk
www.lovibond.com
UK

Tintometer South East Asia

Unit B-3-12, BBT One Boulevard,
Lebu Nilam 2, Bandar Bukit Tinggi,
Klang, 41200, Selangor D.E
Tel.: +60 (0)3 3325 2285/6
Fax: +60 (0)3 3325 2287
lovibond.asia@tintometer.com
www.lovibond.com
Malaysia

Tintometer India Pvt. Ltd.

Door No: 7-2-C-14, 2nd, 3rd & 4th Floor
Sanathnagar Industrial Estate,
Hyderabad, 500018
Telangana
Tel: +91 (0) 40 23883300
Toll Free: 1 800 599 3891/ 3892
indiaoffice@tintometer.com
www.lovibondwater.in
India

Tintometer AG

Hauptstraße 2
5212 Hausen AG
Tel.: +41 (0)56/4422829
Fax: +41 (0)56/4424121
info@tintometer.ch
www.tintometer.ch
Switzerland

Tintometer Brazil

Caixa Postal: 271
CEP: 13201-970
Jundiaí – SP
Tel.: +55 (11) 3230-6410
sales@tintometer.com.br
www.lovibond.com.br
Brazil

Tintometer Spain

Postbox: 24047
08080 Barcelona
Tel.: +34 661 606 770
sales@tintometer.es
www.lovibond.com
Spain